



Nozzle Cleaning Procedure
MAGIC 200
Steven Spielman
Aerosol Dynamics Inc.
2019-06-13

- Obtain a soft interdental brush. Finer size is better. For example, is version by Gum® works:

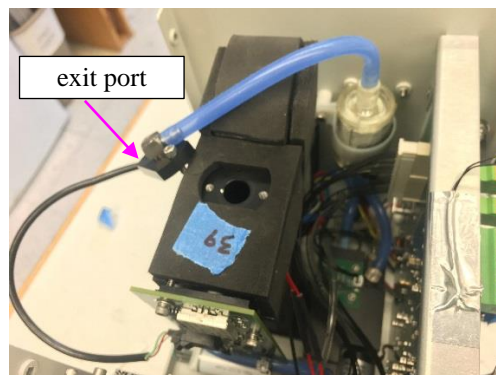

- Power off the unit.
- Remove the lid from the case (8 Phillips-head screws).
The foam-covered part near the top of the interior is the optics head.
- Use a 3/32 hex wrench to remove the two screws securing the exit port at the top of the optics head. Move the exit port aside, being careful not to lose the screws and o-ring.

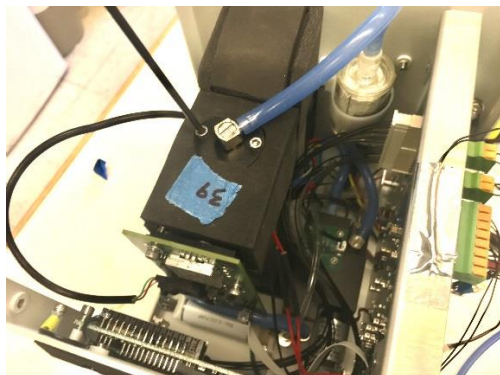

- With the port removed, find the end of the nozzle. It should be visible about 14 mm below the opening, pointing upward.

- Insert the tip of the dental pick into the end of the nozzle. Slide it up and down a couple of times and pull it out.



- Replace the exit port. Make sure the o-ring is in the groove on the underside.
- Attach the case lid.